

FACILITATED AND RELIABLE PROCESS MONITORING DURING VACCINE PROCESS DEVELOPMENT AND PRODUCTION USING THE AUTOMATED SAMPLING SYSTEM NUMERA®

Pascal Vonlanthen Securecell AG, In der Luberzen 29, 8902 Urdorf, Switzerland (pascal.vonlanthen@securecell.ch)

Simon Flückiger Malcisbo AG, Wagistrasse 27a, 8952 Schlieren, Switzerland (info@malcisbo.com)

Lukas Neutsch ZHAW, Grüentalstrasse 14, 8820 Wädenswil, Switzerland (lukas.neutsch@zhaw.ch)

ABSTRACT

Vaccines are one of the most effective, cost-efficient, and sustainable ways of controlling infectious diseases and have had a significant impact on public health and longevity as well as intensive livestock production. Vaccines are generally classified as live or non-live vaccines. Live vaccines contain attenuated replicating strains of pathogens, whereas non-live vaccines contain inactivated whole pathogens or components of pathogens. Over the past years, so-called third-generation vaccines have been developed including RNA and DNA vaccines, viral vector vaccines, and virus-like particles.

In a collaboration between Zurich University of Applied Science and Malcisbo AG, an orally administered, live-attenuated vaccine against *Campylobacter* infections for chickens is being developed, using Malcisbo's glycoengineering platform technology. The highly conserved protein-bound N-glycans of *Campylobacter jejuni* and *Campylobacter coli* are recombinantly expressed on the outer membrane of an attenuated *Salmonella* strain. During a continuous fermentation, different process conditions were tested to optimize antigen expression and cell size distribution of the recombinant *Salmonella* strain. With Securecell's automated sampling system Numera®, the bioprocess was sampled over 245 h, complementing manual sampling and enabling the reliable high-frequency monitoring of critical material attributes.

INTRODUCTION

The use of vaccines has transformed public health and many of the diseases that were previously responsible for the majority of childhood deaths have largely disappeared including Diphtheria, Measles, Pertussis, and Polio [1]. Vaccines are biological products that exploit the ability of the highly evolved mammalian immune system to respond to and remember encounters with pathogens. To induce protective immunity, the vaccine must contain pathogen-derived antigens. Vaccines are generally

classified as live(-attenuated) and non-live vaccines depending on the nature of the antigens. Live-attenuated vaccines contain replicating strains of the relevant pathogen while non-live vaccines contain non-replicating pathogens or components of the pathogen. Vaccines containing components of a pathogen are known as subunit vaccines encompassing protein vaccines, polysaccharide vaccines, or conjugate vaccines. In addition to the traditional live and non-live vaccines, other types of

vaccines have been developed for example nucleic acid-based RNA and DNA vaccines, viral vectors vaccines, and virus-like particles [2]. In most cases, live vaccines elicit a wider range of immunologic responses activating both humoral (B cells) and cellular (CD4+ and CD8+ T cells) reactions than non-live vaccines and therefore also have significantly higher immunogenicity. Often, a single vaccine administration is sufficient to induce long-term protection [3].

Vaccines have not only transformed public but also veterinary health and are widely accepted as a cost-effective and sustainable immunization method for companion animals and livestock to efficiently prevent the transmission of zoonotic diseases [4]. *Campylobacter* is one of the most important foodborne pathogens in Europe and the US, and the causative agent of campylobacteriosis. The disease manifests itself through mild to severe inflammatory diarrhea and can result in chronic implications such as reactive arthritis or Guillain-Barré-Syndrome. The main transmission occurs through the consumption of contaminated poultry meat which causes about 60 - 80% of global campylobacteriosis cases. Frequently, isolated *Campylobacter* from human infections showed resistance to various antibiotics. There are at present no vaccines approved by any global governing authority to prevent *Campylobacter* infections. Thus, vaccines represent a major unmet medical and industrial need to reduce the use of antibiotics, prevent the spread of antibiotic resistance, and raise food security [5,6].

Zurich University of Applied Science and Malscibo AG are currently developing an orally administered,

live-attenuated vaccine against *Campylobacter* infections for broiler chickens. The high variability of *Campylobacter jejuni* (main *Campylobacter* gastroenteritis-causing member) with more than 60 serotypes makes vaccine generation against all serotypes a challenging endeavor [7]. Using Malscibo's glycoengineering platform technology a vaccine candidate covering all serotypes was established. The highly conserved protein-bound N-glycans of *Campylobacter jejuni* and *Campylobacter coli* are recombinantly expressed on the outer membrane of an attenuated *Salmonella* strain. First experiments using the glycoengineered vaccine against *Campylobacter* in chicken showed a reduction of *Campylobacter jejuni* in the chicken feces, suggesting a good efficacy of the vaccine candidate [8].

Process development and optimization

Over the last decade, cost control in biopharmaceutical manufacturing has become more relevant as margins, especially for vaccines compared to other biopharmaceutical products, are low. It is important to decide early in the development process how the product will be manufactured at a commercial scale. Changes later on such as new manufacturing equipment or raw material components will typically trigger new regulatory requirements, including clinical trials and thus additional costs [9].

To make an early decision on the process details i.e., optimal process conditions for the manufacturing of the *Campylobacter* vaccine candidate, the substrate concentration of critical growth medium components was altered marking different process phases

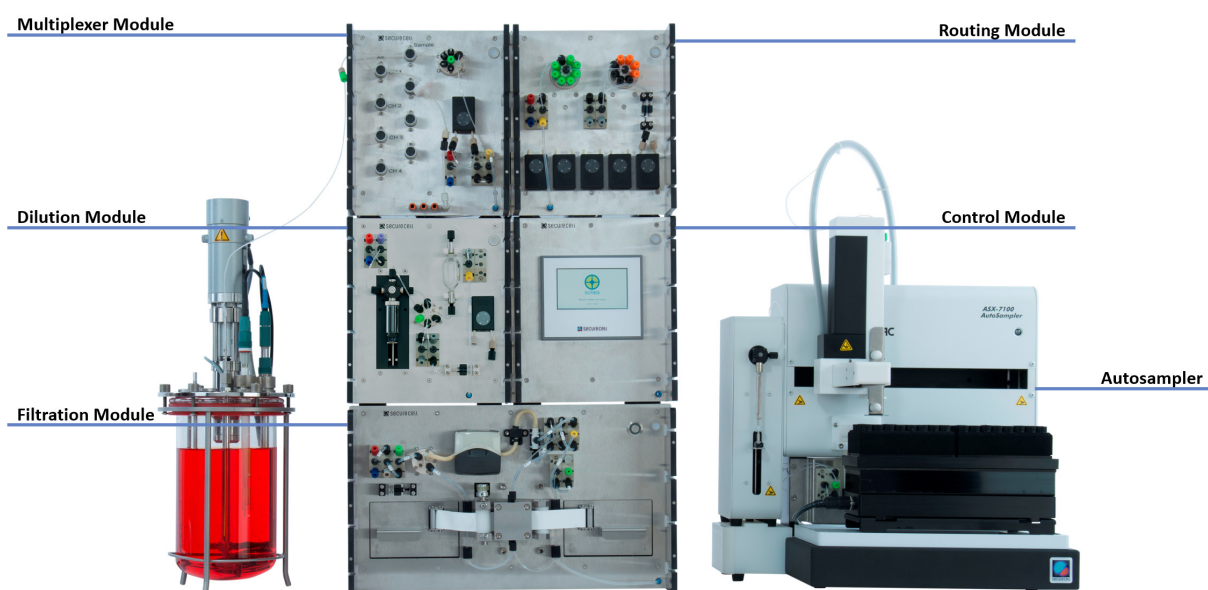


Figure 1: Image of the Numera® automated sampling system. The Multiplexer Module forms the interface to the bioreactors. After optional sample processing by the Dilution or Filtration Module, the Routing Module directs the sample either to the autosampler for sample storage and off-line analysis or directly to connected 3rd party analyzers.

during continuous culture and the influence on the antigen expression (AE) and cell size distribution was evaluated. To be able to track and reliably measure the surface AE and cell size distribution of the *Salmonella* culture with flow cytometry, liquid samples were taken in regular intervals for 245 h from the bioreactor and were analyzed. Securecell AG's modular sampling system Numera® was used to complement manual sampling outside working hours by automatically drawing and storing samples in a cooled environment, making samples readily available for flow cytometric analysis.

Numera® can be flexibly configured to allow sampling of up to 16 bioreactors or downstream devices in parallel, seamless in-built sample processing, and sample transfer to a sample storage unit or 3rd party analyzers (Figure 1). Compared to other sampling systems, Numera® allows for sampling of small sample volumes at a high frequency, reliable sample processing with a precise dilution (error $\leq 2\%$ SD) and unique tape-filtration technology, cooled sample storage, and well-established integration with multiple cell and biochemical analyzers and HPLC systems.

Material and Methods

All experiments described were performed in a bioprocessing laboratory with a fully integrated IoT infrastructure (i²BPLab) at Zurich University of Applied Sciences (ZHAW), Wädenswil.

Cultivation conditions

The total process duration was 245 h including a 6h preparation phase for a sterility check and media conditioning phase.

The continuous fermentation was performed in a 1.5L Minifors 2 bioreactor (Infors HT) equipped with a pH

and a dissolved oxygen (DO) sensor. Additionally, the reactor was equipped with a feed tube and two dip tubes. One dip tube was connected to an external peristaltic pump to set the culture volume during the continuous phase (F_{out}). The second dip tube was split by a Y-connector and connected to Super Safe Samples (Infors HT), for manual sampling, and directly to the Numera® Multiplexer Module, for automated sampling.

The reactor was supplied with air (1.5 L min⁻¹) and O₂ to control dissolved O₂ around minimally 25 %. The temperature and stirrer setpoints were set to 37 °C and 1200 rpm, respectively. The pH was controlled at 7.2 by the addition of 25 % NH₄⁺ and 8.5 % H₃PO₄.

Set-up automated sampling

The reactor system and the Numera® system were connected to and controlled by the software Lucillus® (Securecell AG). In Lucillus®, the sampling events were planned and, in the case of automated Numera® sampling, triggered during the process execution phase. From the samples, the surface AE level, the cell size distribution of the population, and the optical density at wavelength 600 nm (OD₆₀₀) were determined. Additionally, for the manually drawn samples only, the cell dry weight (CDW) and the substrate concentration in the culture medium were measured. Other process-related parameters were measured in-line with the respective probes. Sampling intervals ranged from 1 – 6 h depending on the process phase.

Off-line sample analysis

CDW was measured by pelleting 2 mL of culture broth in a pre-weighed tube, discarding the supernatant, and drying the pellet at 105 °C for 48 h. For the estimation of the AE and cell size, the cells were

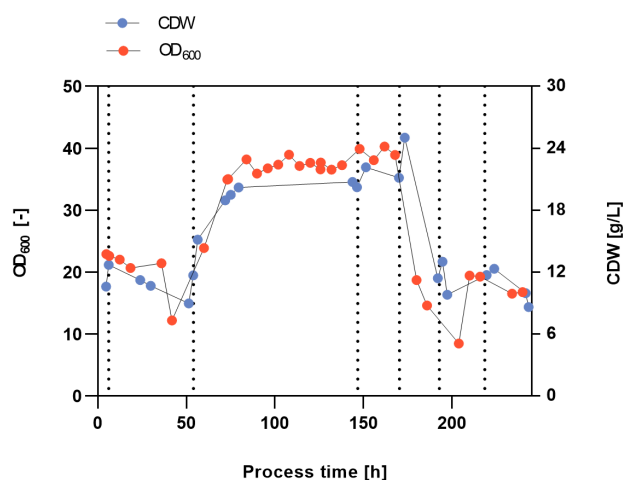


Figure 2: CDW measured after manual reactor sampling and OD600 measured after automated Numera® sampling over the six different process phases indicated by vertical dashed lines with varying substrate concentrations.

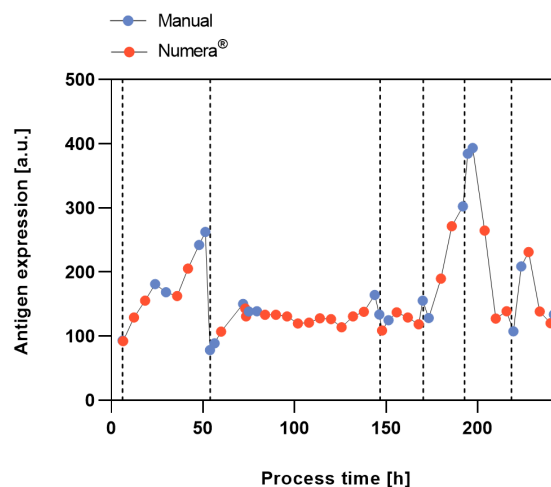


Figure 3: Antigen expression after manual or automated Numera® sampling over the six different process phases indicated by vertical dashed lines with varying substrate concentrations.

stained with a fluorescent stain and subsequently analyzed by flow cytometry.

Results and Discussion

To explore optimal process conditions for the manufacturing of the vaccine candidate, the glycerol, tryptone, and yeast extract ratio in the growth medium was altered marking different process phases indicated by vertical dashed lines during the continuous culture (Figure 2-4). The aim of Malsisbo's scientists was on the one hand process development and optimization (i.e., cell density) for upcoming industrial *Campylobacter* vaccine production, and on the other hand vaccine quality optimization (i.e., maximization of AE and cell size). Larger cells have a larger surface area and can thus express more of the antigen.

The cell densities over the complete process were measured after manual sampling and heating (CDW) or after automated sampling with the Numera® sampling system (OD_{600}) (Figure 2). The cell density varied considerably in the different process phases most likely due to the different substrate concentrations. During phase 1, an unintended washout of the cells occurred from which the culture recovered. Both manual and Numera® automated cell density measurements showed a similar growth curve over the complete process. The cell density peaked at the start of phase 4.

The manual and Numera® samples were also analyzed by flow cytometry to determine the surface AE (Figure 3) and relative cell size distribution of the population (Figure 4) over time. Also, the level of AE

altered considerably in the different process phases. The increase of AE during process phase 4 was most pronounced, followed by phase 1.

Similarly, the relative cell size distribution (FSCI-1 – FSCI-10) changed over process time. The relative number of big cells (> FSCI-6) was highest in mid-phase 4, followed by mid-phase 1.

Conclusion

Understanding how the cell density, the antigen presentation, and the relative cell size distribution are influenced by the process conditions is crucial.

From a supplier perspective, Numera® successfully complemented manual sampling. An unattended Numera® sampling run between 80 h and 145 h process time delivered consistent results for OD, AE, and cell size. This exemplarily demonstrated the reliability of the Numera® system for high-frequency sampling during non-office hours and storage of the samples for later (flow cytometric) analysis.

From a scientific perspective, these preliminary and exploratory experiments showed that substrate concentration for AE and cell size seemed to be optimal during phases 1 and 4 as in these phases AE and the relative number of big cells (> FSCI-6) peaked. However, the highest cell densities were reached during phase 2. This observation might be explained by the fact that after the washout, the cells invested energy and resources in proliferation instead of surface antigen expression. As the influence of the washout cannot be fully accounted for, the results have to be verified in subsequent experiments.

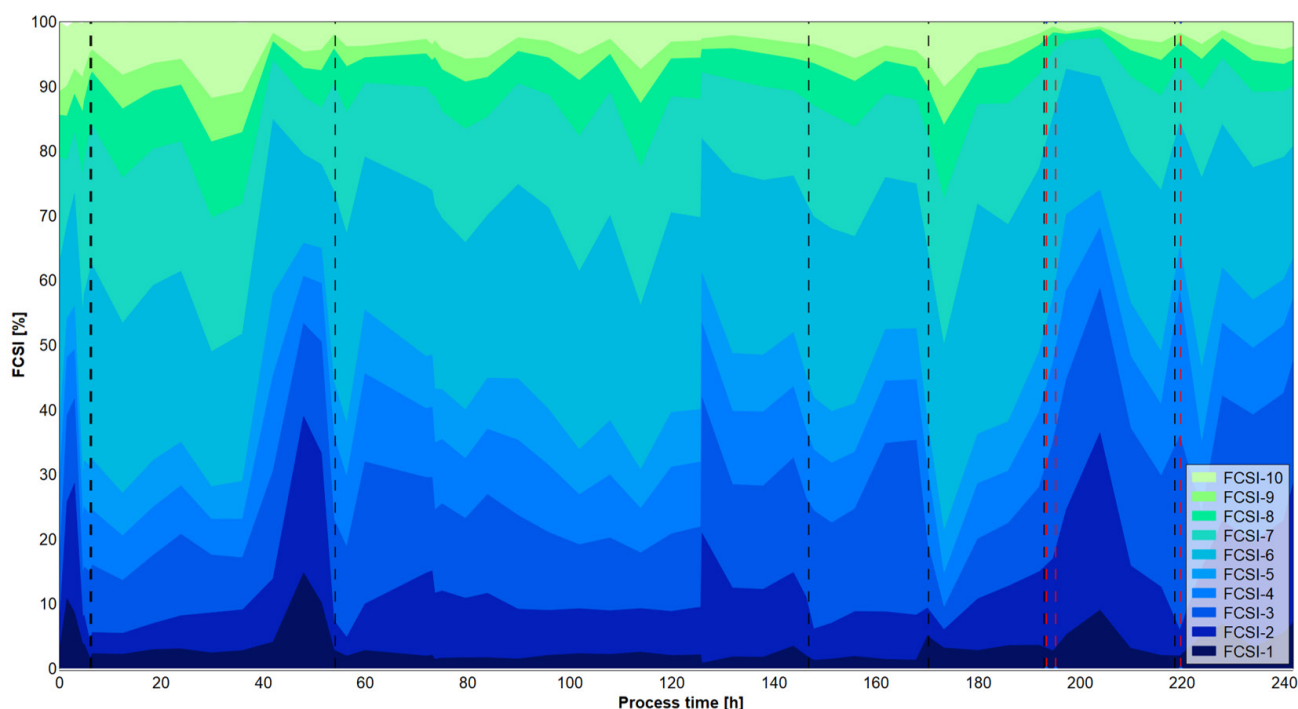


Figure 4: Cell size distribution measured after automated Numera® sampling over the six different process phases indicated by vertical dashed lines with varying substrate concentrations. FSCI-1: Very small cells. FSCI-10: Very large cells.

KEY POINTS

- Highly frequent and automated bioreactor sampling and sample storage with Numera®
- Reliable sampling during non-office hours
- Good accordance of off-line and on-line analytical results
- Combination of Numera® and Lucillus® as integrated PAT solutions
- Complete automation of sample-based analytics

REFERENCES

1. World Health Organization. Global vaccine action plan 2011-2020. WHO https://www.who.int/immunization/global_vaccine_action_plan/GVAP_doc_2011_2020/en/ (2013)
2. Pollard, A. J. & Bijker, E. M. A guide to vaccinology: from basic principles to new developments. Nat Rev Immunol 21, 83-100 (2021)
3. Kollaritsch, H. & Rendi-Wagner, P. 9 - Principles of Immunization. in Travel Medicine (Third Edition) (eds. Keystone, J. S., Freedman, D. O., Kozarsky, P. E., Connor, B. A. & Nothdurft, H. D.) 67-76 (Elsevier, 2013)
4. Jorge, S. & Dellagostin, O. A. The development of veterinary vaccines: a review of traditional methods and modern biotechnology approaches. Biotechnology Research and Innovation 1, 6-13 (2017)
5. Igwaran, A. & Okoh, A. I. Human campylobacteriosis: A public health concern of global importance. Heliyon 5 (2019)
6. Facciola, A. et al. Campylobacter: from microbiology to prevention. J Prev Med Hyg 58, E79-E92 (2017)
7. Munroe, D. L., Prescott, J. F. & Penner, J. L. Campylobacter jejuni and Campylobacter coli serotypes isolated from chickens, cattle, and pigs. J Clin Microbiol 18, 877-881 (1983)
8. <https://www.malcisbo.com/technologies-projects#campylobacter> (2022)
9. Plotkin, S., Robinson, J. M., Cunningham, G., Iqbal, R. & Larsen, S. The complexity and cost of vaccine manufacturing - An overview. Vaccine 35, 4064-4071 (2017)

© Copyright 2022 by Securecell AG. All rights reserved, including graphs and images. Lucillus and Numera are registered trademarks of Securecell AG, Switzerland



SECURECELL AG
In der Luberzen 29
8902 Urdorf, Switzerland

+41 44 732 90 70
contact@securecell.ch

For more information on
our products and services
visit our website at
www.securecell.ch

